



## Flash - Dos and Don'ts

### Do

- Clean arc pins, sample sensor and oven surface
- Keep your sample cups free of dings or scratches
- Schedule recommended service/PM
- Perform daily or as needed "Bakeout" - "Daily Maintenance Procedure"

#### 4. DAILY MAINTENANCE

---

To guarantee repeatable results the daily maintenance procedure should be performed.


Please clean the surface of the oven plate well with tissue and solvent. do not **scratch**.

The surface must be free of residuals and scratches for good heat contact.


Do not bend the temperature sensor and the arc pins.

Clean the arc pins (1) and the temperature sensor (2) carefully with tissue and the brass eraser.

Cleaning the arc pins is necessary to regain a metal surface at the tips for easy ignition.

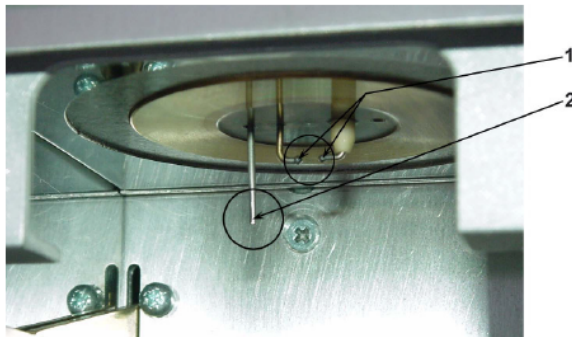

FL-RADIERER-MESSING \ Brass eraser

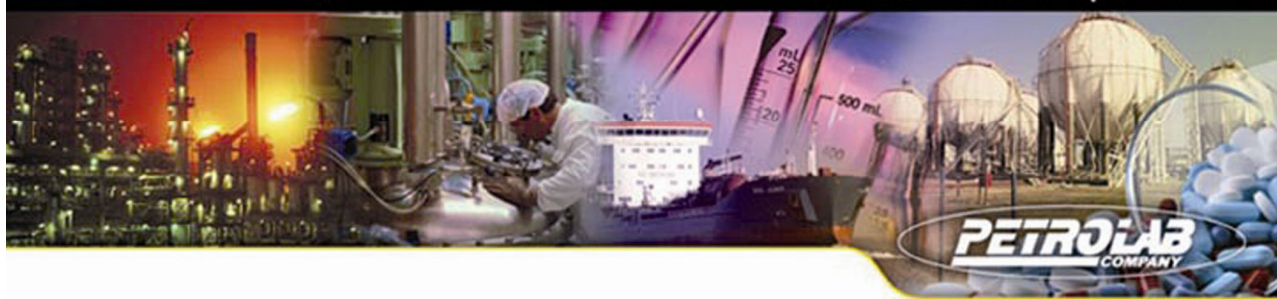

### Don't

- Let sample buildup on oven surface or arc pins
- Throw sample cups on top of each other
- Neglect regular service intervals - Factory Recommendations
- Bend or damage sample sensor or arc pins

### Don't Examples

